

IMPORTANT INFORMATION – PLEASE READ

Installation Requirements for WASP Touch

Thank you for choosing one of our products for your laboratory. To enable our engineers to perform an efficient, trouble-free installation please study, complete and return this form to us at: service@dwscientific.co.uk.

Should you have any questions, please do not hesitate to contact us, as we are here to help. When we have received the completed form, our Service Department will contact you to arrange a mutually convenient installation date.

Please contact us if you have any further questions about the installation procedure.

Space Requirements

The weight of the equipment is 21kg. The bench allocated must be flat, level and of sufficient size to support the base fully.

External Dimensions: WASP Touch

1 litre Sanitizing Station

2 litre Sanitizing Station

Width	Depth	Height		Width	Depth	Height		Width	Depth	Height
<i>mm</i>	<i>mm</i>	<i>mm</i>		<i>mm</i>	<i>mm</i>	<i>mm</i>		<i>mm</i>	<i>mm</i>	<i>mm</i>
420	374	298		120	330	270		152	412	320

1. Have you allocated flat, level bench space that can support 21kg and an item of 420 x 374 x 298mm (D x L x H) and the Sanitizing System and power supply?
2. Have you allocated a spare mains outlet for the unit (230 +/- 10%V AC)?
3. Do you have suitable disinfectant available to fill the bottles?
4. Have you seen Technical Note WA06 which sets out the best way to prepare agar plates for the WASP?
5. Have you ordered sufficient disposable beakers?
6. Are you fully aware of the Quality Counts Scheme and what actions are required of you (UK and Europe only)? Yes / No
7. Will real samples be put through the WASP on the day of installation? Yes / No
8. If you are manually counting plates, is any additional training required? Yes / No
9. Are you using pre-poured media? Yes / No
If so, who is the supplier?

In the UK, delivery and installation are free of charge (unless otherwise agreed). If our engineers are unable to install the unit and a return journey is necessary, **a charge may be made.**
Export customers should please refer to their local distributor.

It is essential that this form is completed and returned, to avoid delay to your installation.

THANK YOU FOR THINKING WHITLEY

Signature

Title

Print Name

Establishment